

Old Separations on New Resins and New Separations on Old Resins

Phil Horwitz and Dan McAlister

PG Research Foundation, Inc.

8205 S. Cass Ave., Suite 106, Darien, IL 60561



Rapid Separations Designed for Automation

Soil 100mg, Water 10mL

Separation Options

Systems

Actinides (Th, U, Np, Pu, Am)

TEVA/TRU

Sr, Tc (Th, U, Pu, Am)

WBEC/UTEVA/Crown-DGA



Sr, Tc (Np, Pu, Am)

WBEC/UTEVA/Crown-DGA

WBEC = Weak Base EC Resin


Weak Base EC Resin

40% alamine-336 (tertiary amine) on inert support

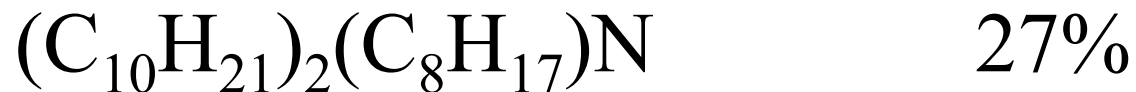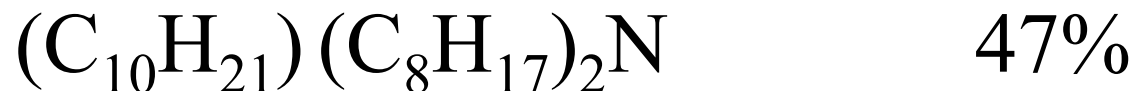


Equilibria for Pertechnetate on Weak Base EC Resin

Extraction:

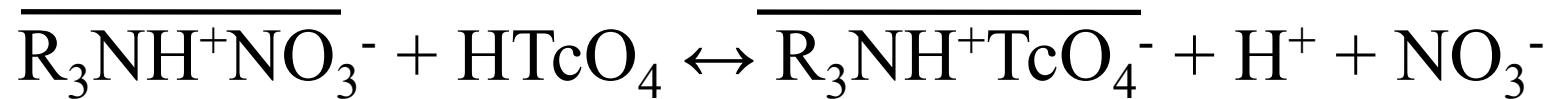


Stripping:

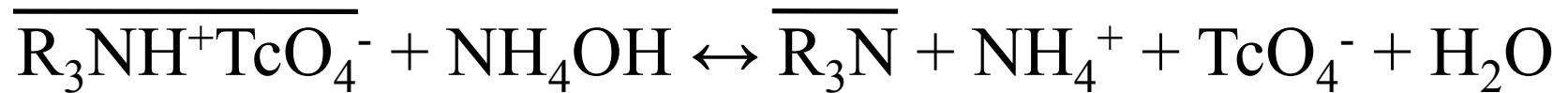


k' on Weak Base EC Resin

100-150 μm , 22(1) $^\circ\text{C}$, 1 hr equilibration

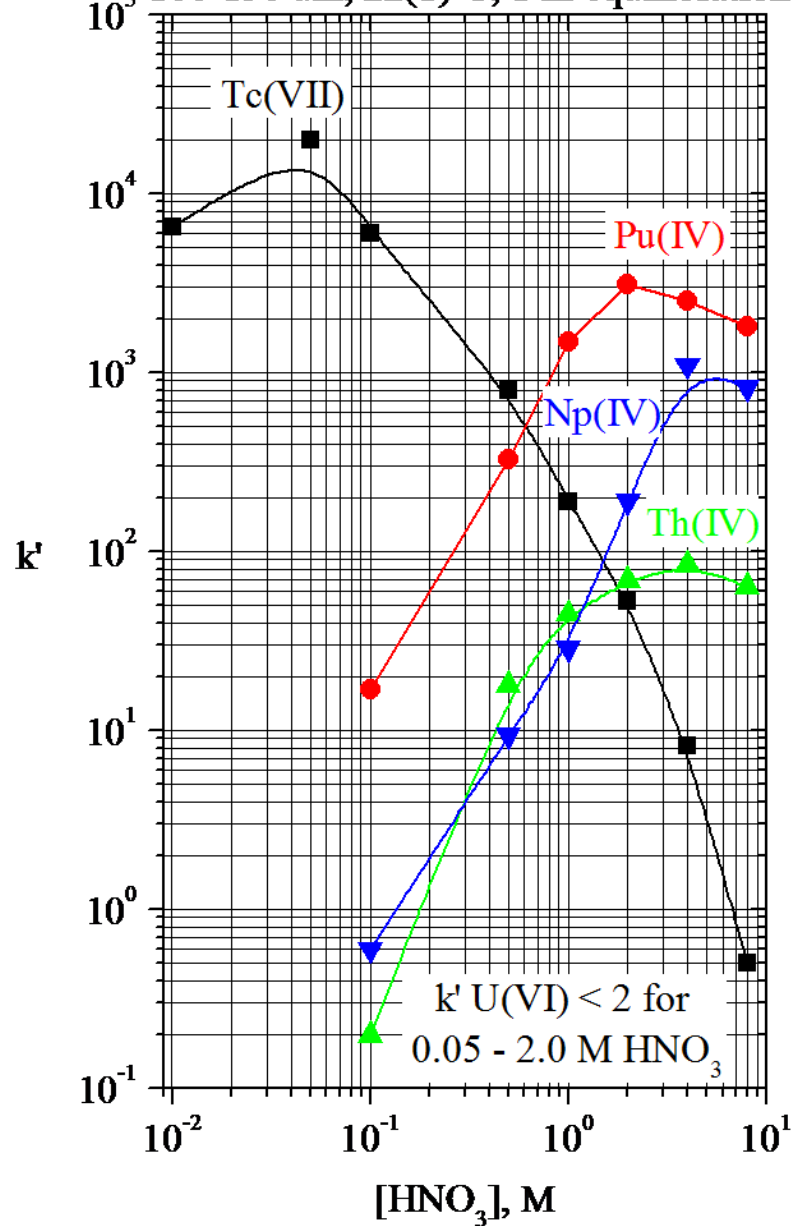


k' on Weak Base EC Resin

50-100 μm , 22(1) $^\circ\text{C}$, 1 hr equilibration

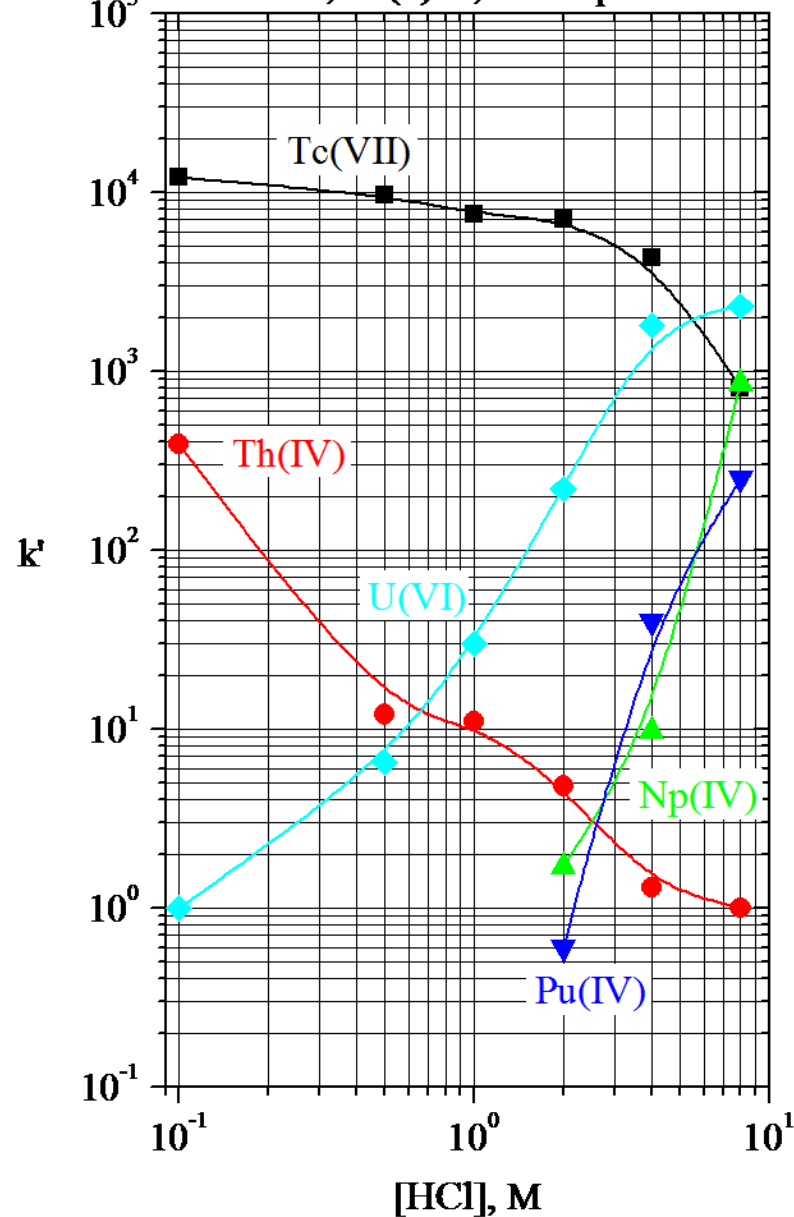


Crown/DGA Resin

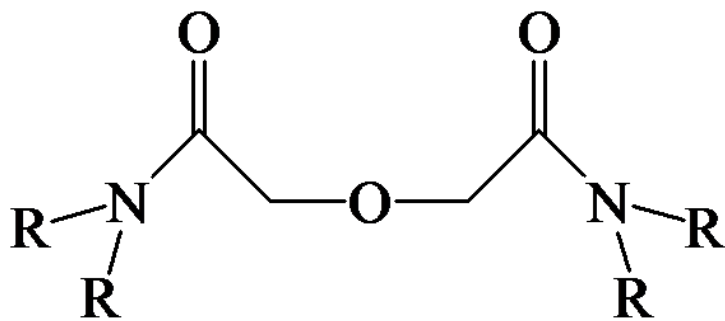

TODGA (R = *n*-octyl)
Ln(III), An(III) (IV) (VI)



DtBCH18-crown-6 present at
the same loading as Sr Resin

TODGA replaces 1-octanol
as diluent for the crown
ether.

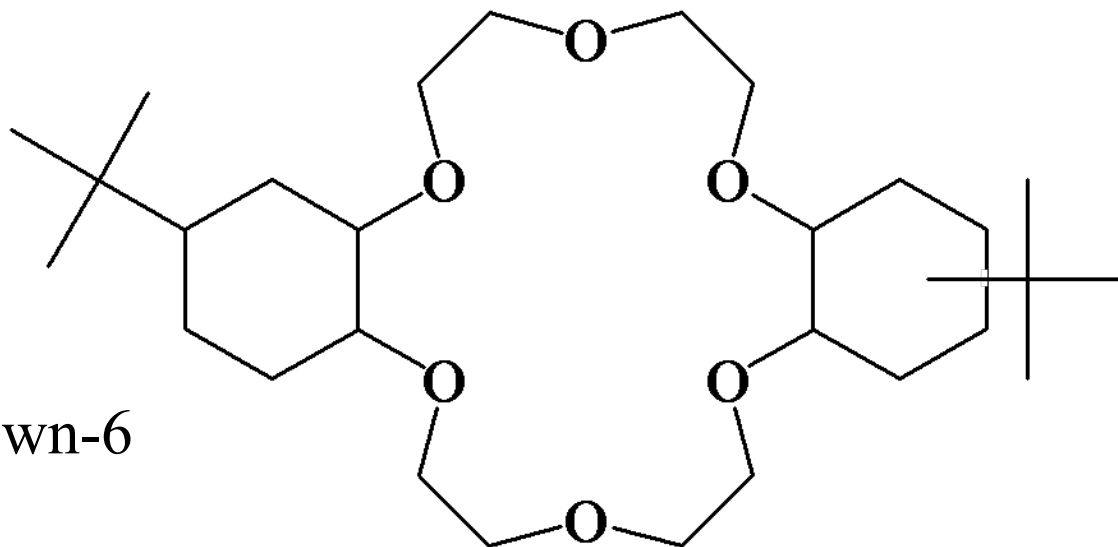


DtBCH18-crown-6
Sr(II)

Equilibria for Extraction of Sr(II) and Am(III)

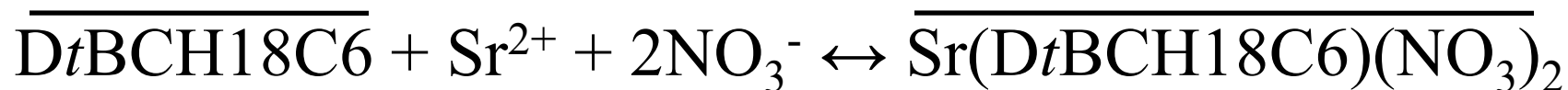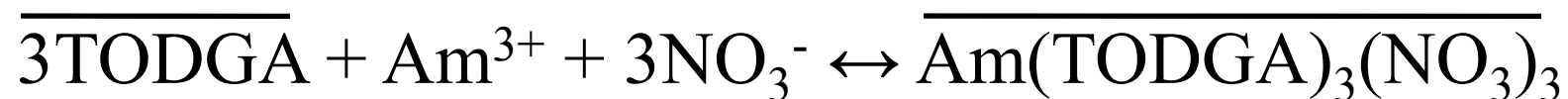


k' Sr(II) on Sr and Crown/DGA Resins

50-100 μm , 22(1) $^\circ\text{C}$, 1 hr equilibration

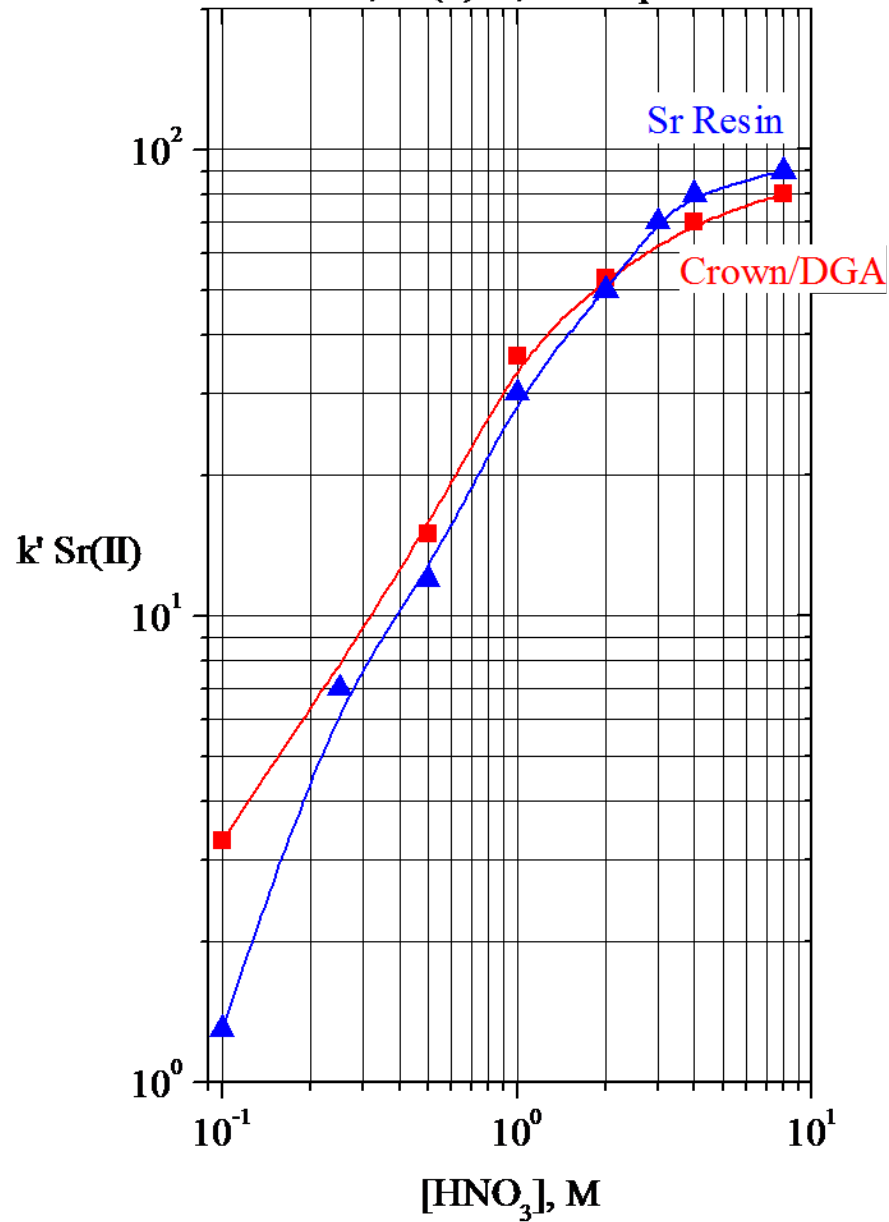


k' Am(III) on DGA and Crown/DGA Resin

50-100 μm , 22(1) $^\circ\text{C}$, 1 hr equilibration

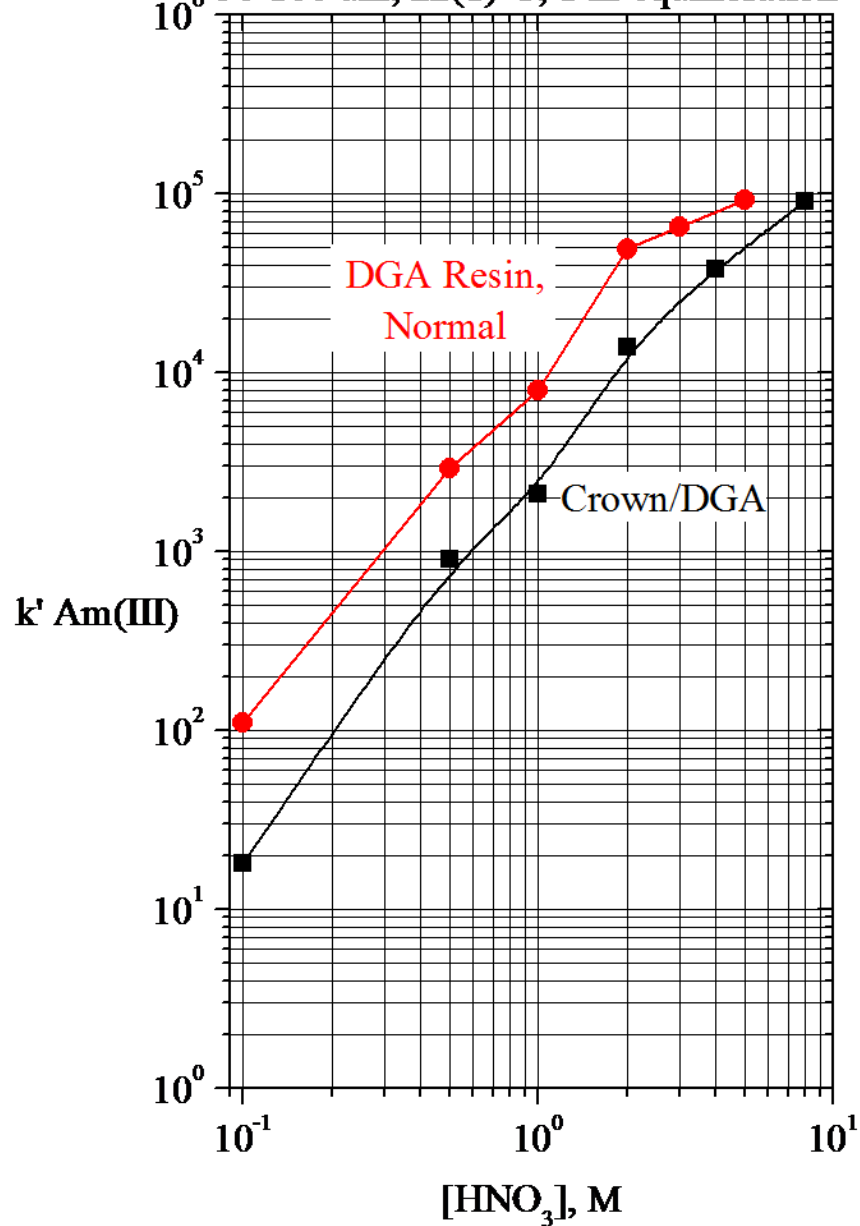


k' on Crown/DGA Resin vs. HNO_3


50-100 mm, 1 h contact time, 22(1) $^\circ\text{C}$

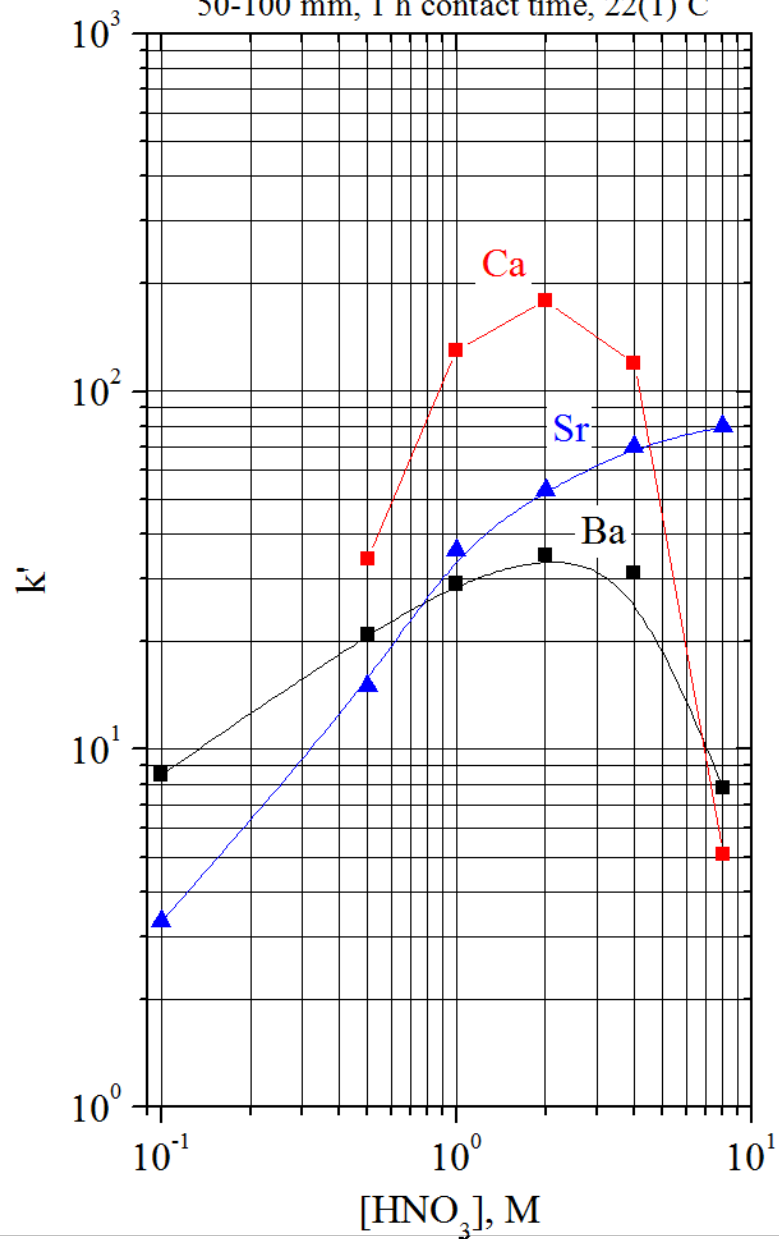

k' Am(III) on DGA and Crown/DGA Resin

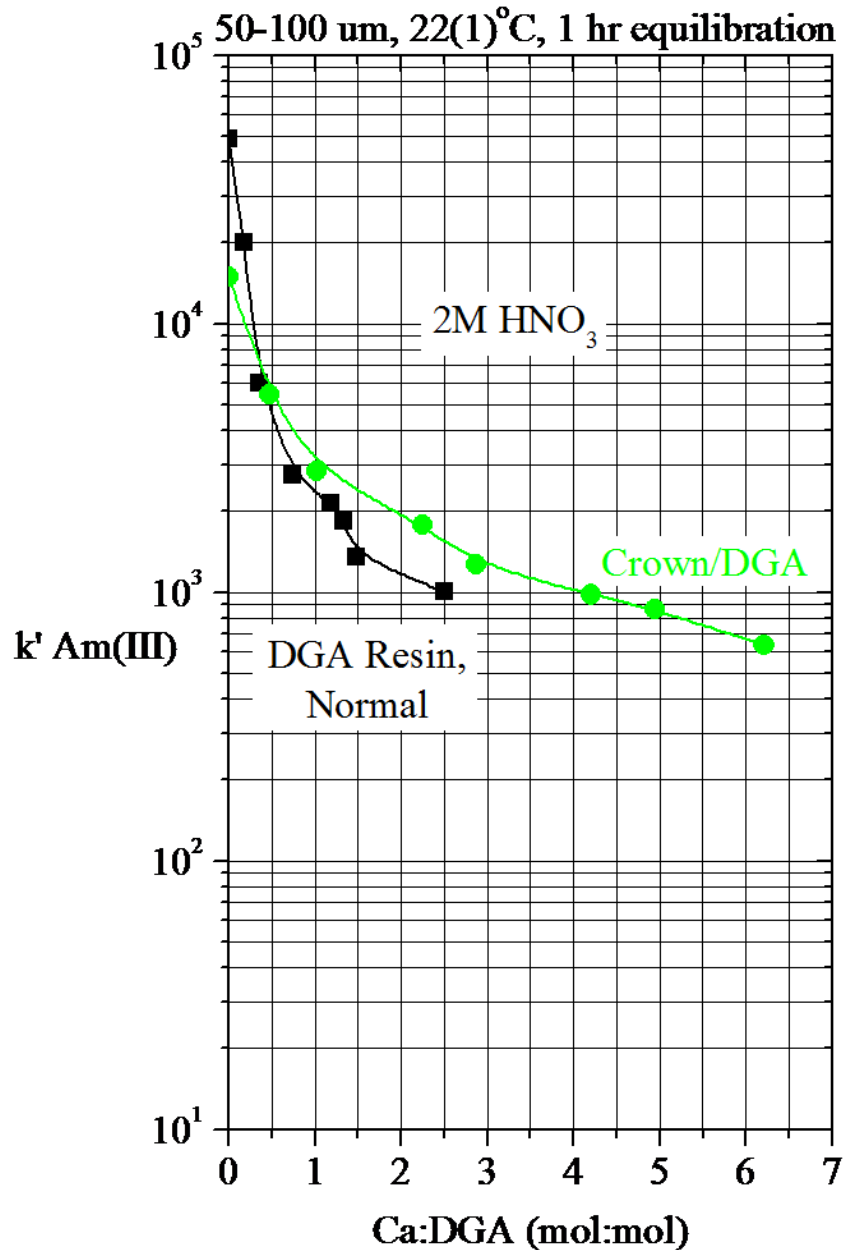

k' Sr(II) on Sr and Crown/DGA Resin

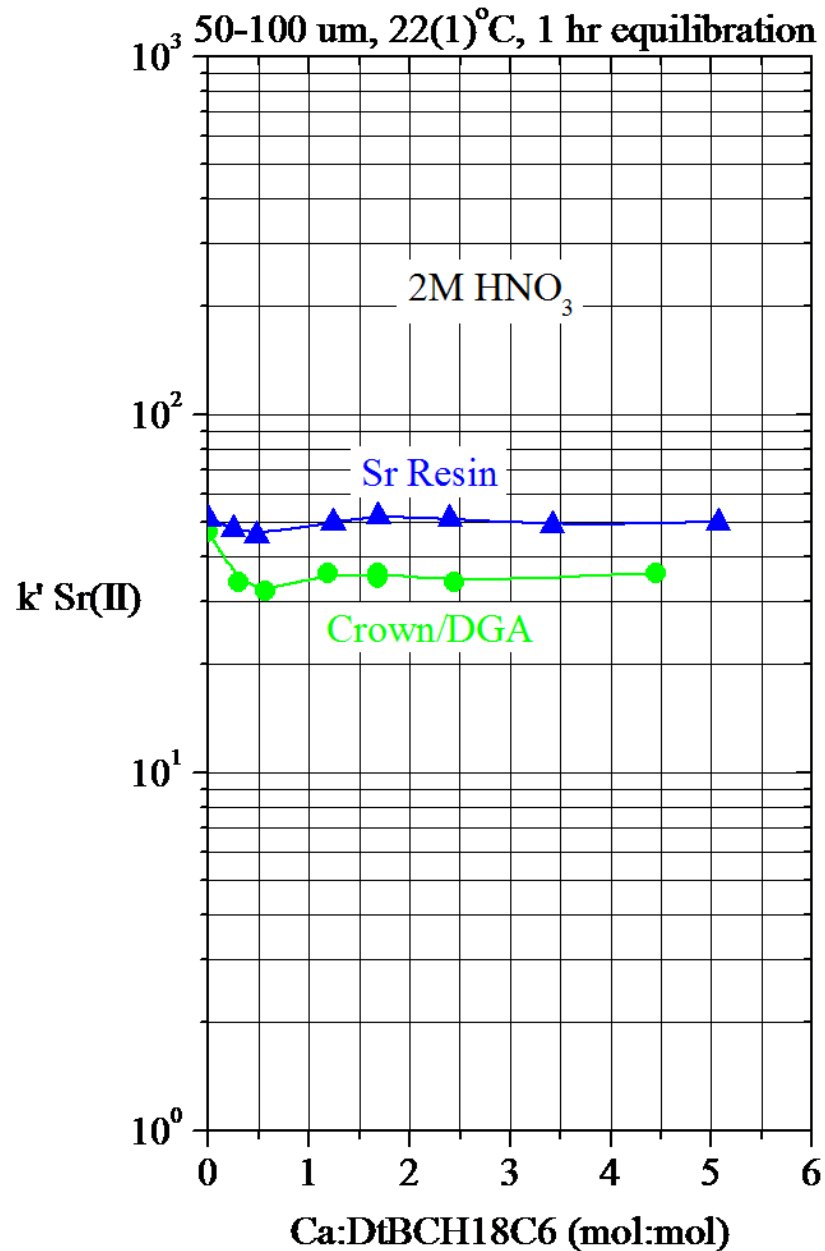

Soil Leachate or Water Sample



acidify to 2M HNO₃

Weak Base EC → Th, Pu, Tc



UTEVA → Th, U



Crown-DGA → Sr, Am





Waste

Separation of Tc, Pu, Th, U, Am and Sr using 2 mL Cartridges (Weak Base EC/UTEVA/Crown-DGA)

Fraction	Solution	BV	Tc	Pu	%			
					U	Th	Sr	Am
Load	2 M HNO ₃ /0.1 M NaNO ₂	5	4.8	0.4	0.2	0.3	0.4	0.2
		5	2.1	0.1	0.1	0.2	0.4	0.2
Rinse	2 M HNO ₃	5	0.0	0.0	0.1	0.3	0.6	0.2
		5	0.0	0.0	0.1	0.3	0.6	0.0
Th Strip	6 M HCl	6	0.2	0.0	0.1	30.1	0.0	0.1
		6	0.2	0.0	0.1	40.0	0.0	0.2
Pu Strip	1 M HCl	6	0.1	98.4	0.3	1.2	0.1	0.6
		6	0.1	99.0	0.2	3.1	0.1	1.6
Tc Strip	1 M NH ₄ OH	6	94.1	0.2	0.3	0.0	0.1	0.0
		6	96.8	0.2	0.3	0.0	0.1	0.0
U Strip	1 M HCl	6	0.4	0.5	98.6	66.7	0.0	0.2
		6	0.5	0.3	98.8	54.1	0.1	0.5
Sr Strip	4 M HCl	6	0.0	0.0	0.2	0.2	97.9	0.0
		6	0.0	0.0	0.2	0.2	97.8	0.0
Am Strip	1 M HCl	6	0.0	0.2	0.1	0.2	0.9	98.4
		6	0.1	0.2	0.1	0.1	0.9	97.3

DGA·FeCl₃ Resin / HCl System

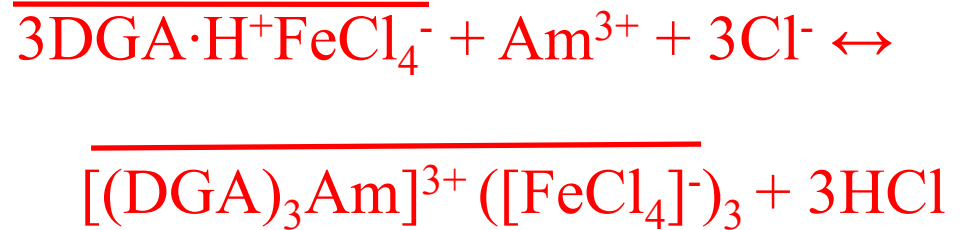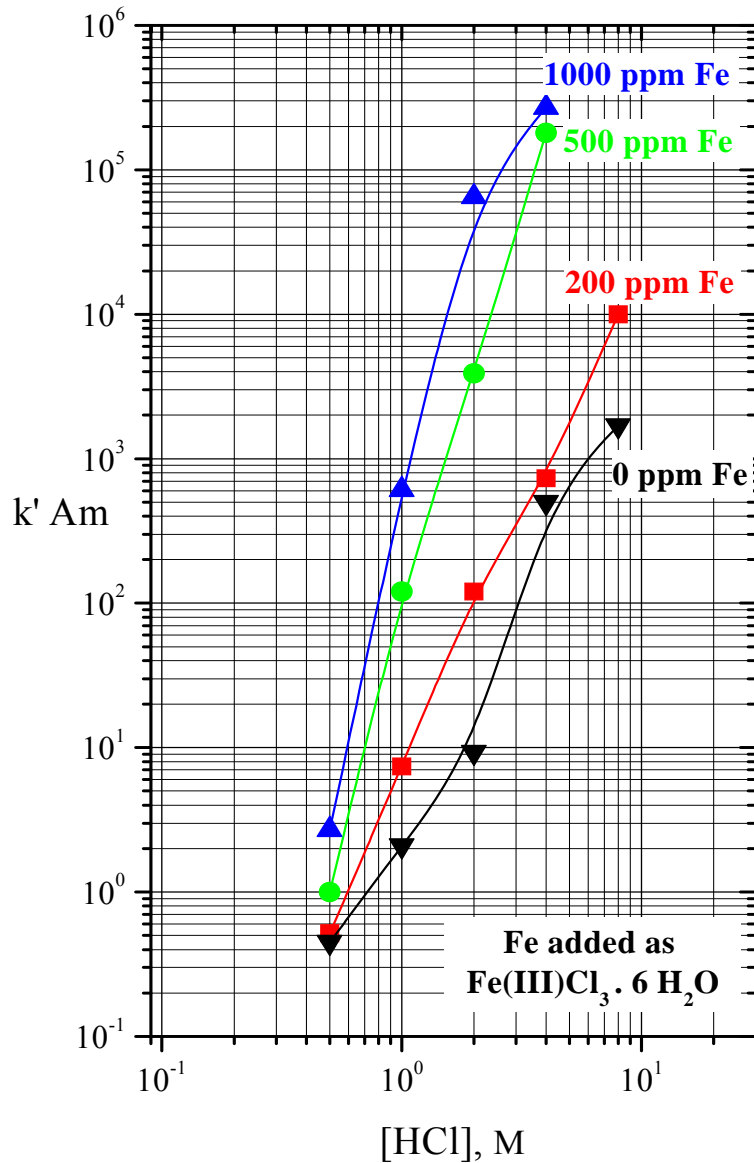

Recovery of Np, Pu and Am from Ferric Hydroxide Precipitate

Ferric Hydroxide Precipitate

↓ 6M HCl

① ~2M HCl Solution + reducing agent

← ② 3M HNO₃


← ③ 0.1M HCl + 0.05M Oxalic Acid

2mL DGA Resin,
Normal Cartridge

Fe + Impurities ← ①②


③ Np, Pu, Am to TRU column


Recovery of ^{227}Ac

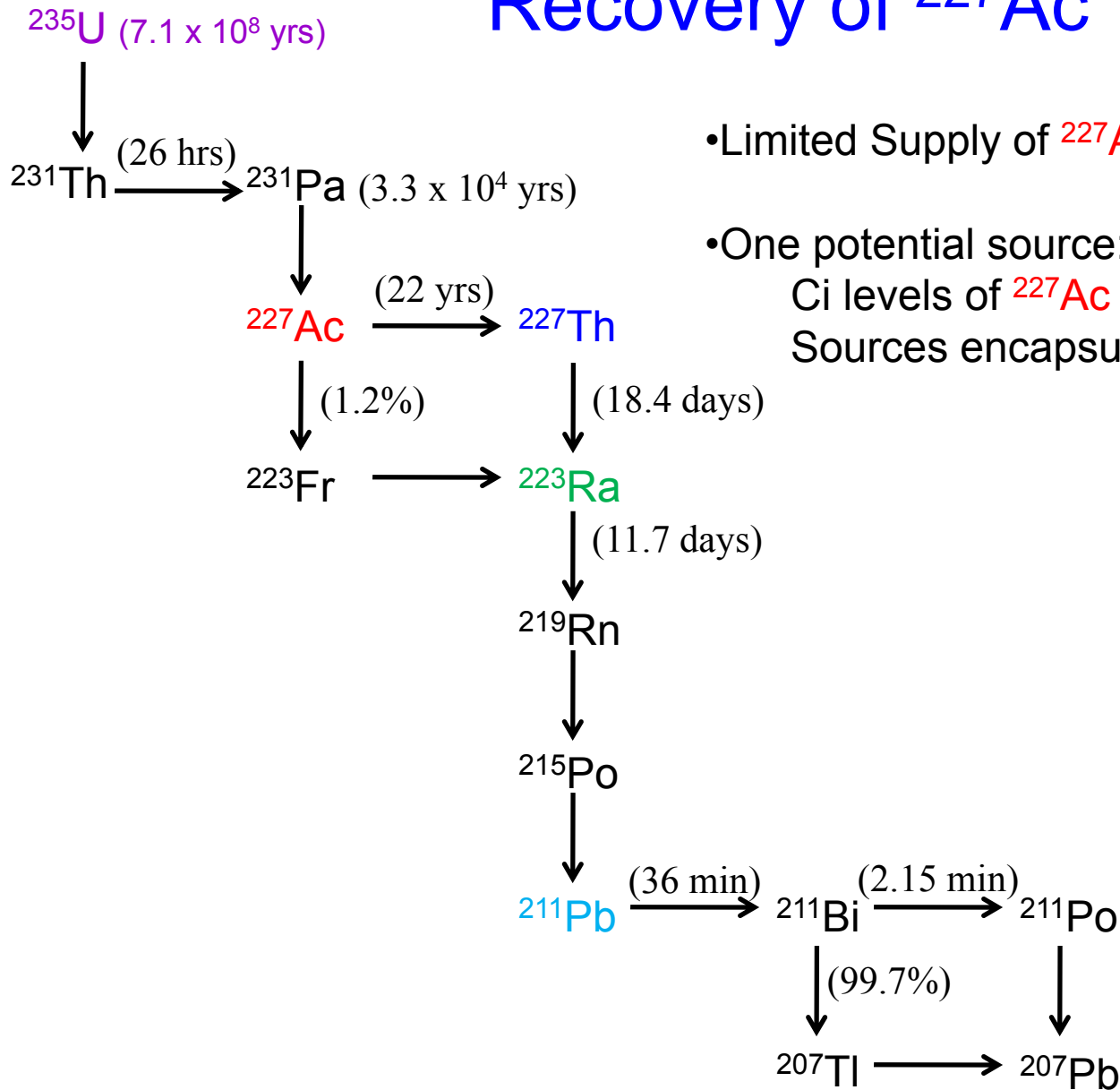

- Limited Supply of ^{227}Ac ($t_{1/2} = 21.77$ yrs)
- One potential source:
Ci levels of ^{227}Ac in Ac/Be Neutron Sources encapsulated with stainless steel

Ac/Th in HCl on 2mL DGA Resin, Normal cartridge

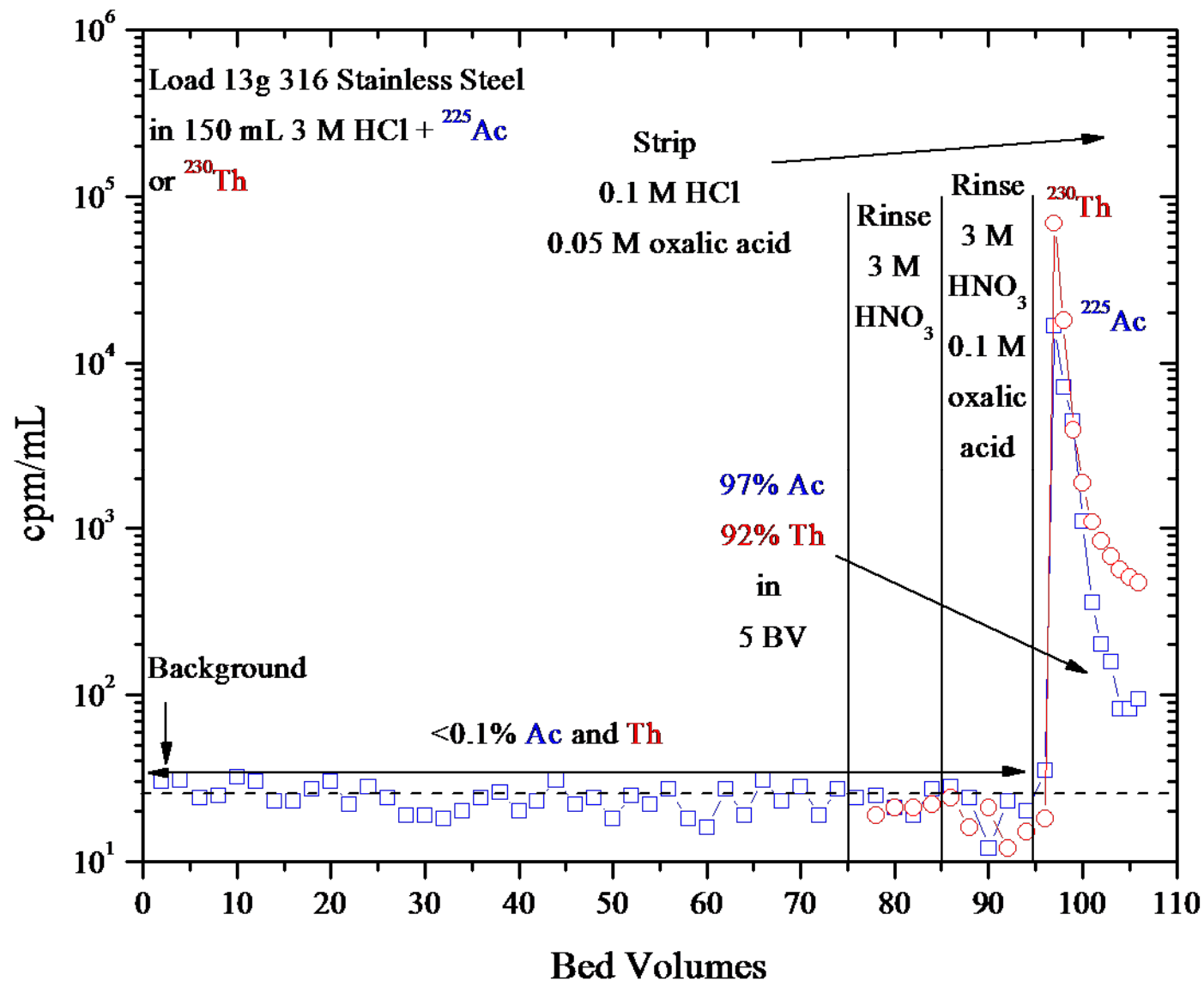

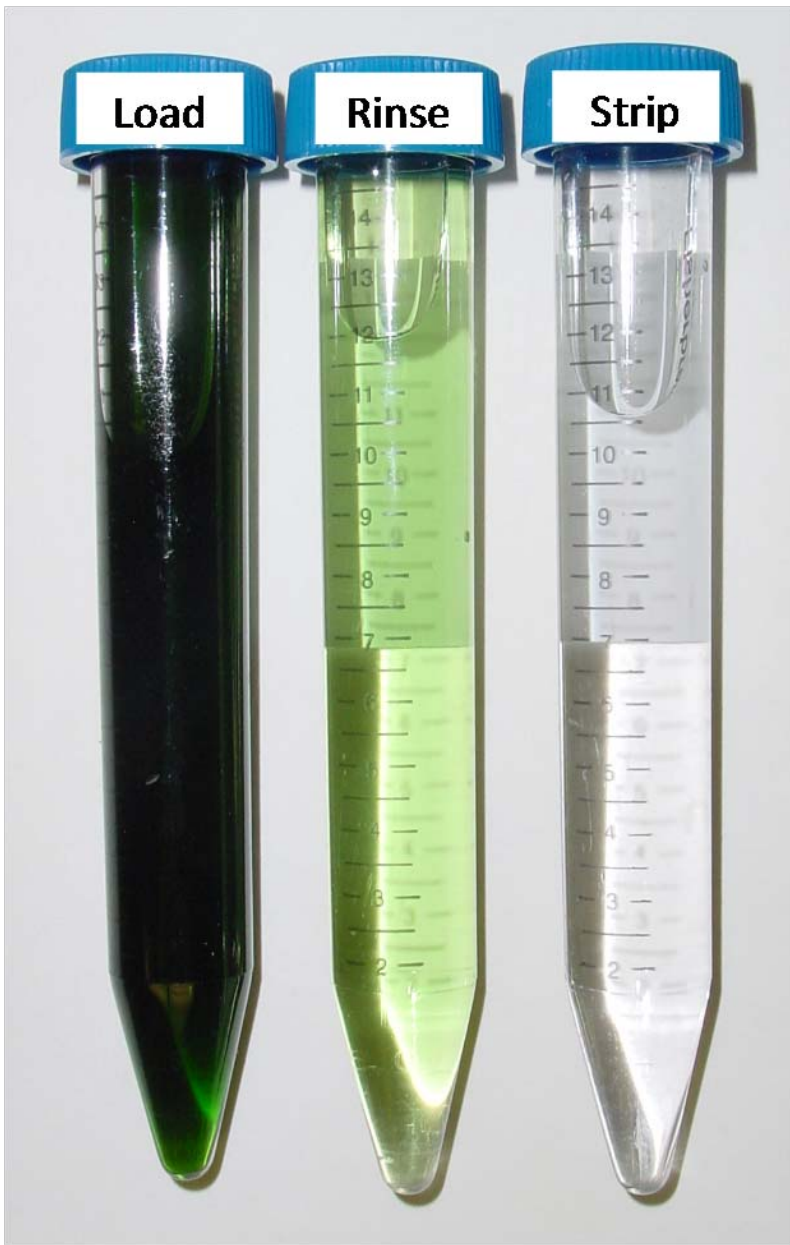

Prefilter

DGA




Recovery of Trace Th from Bulk U

- Applications:

- ^{229}Th (1.4 kg of ^{233}U contains 171 mg (37 mCi) of ^{229}Th)

- ^{234}Th (50 g of ^{238}U contains 17 μCi of ^{234}Th)

Recovery of Trace Th from Bulk U

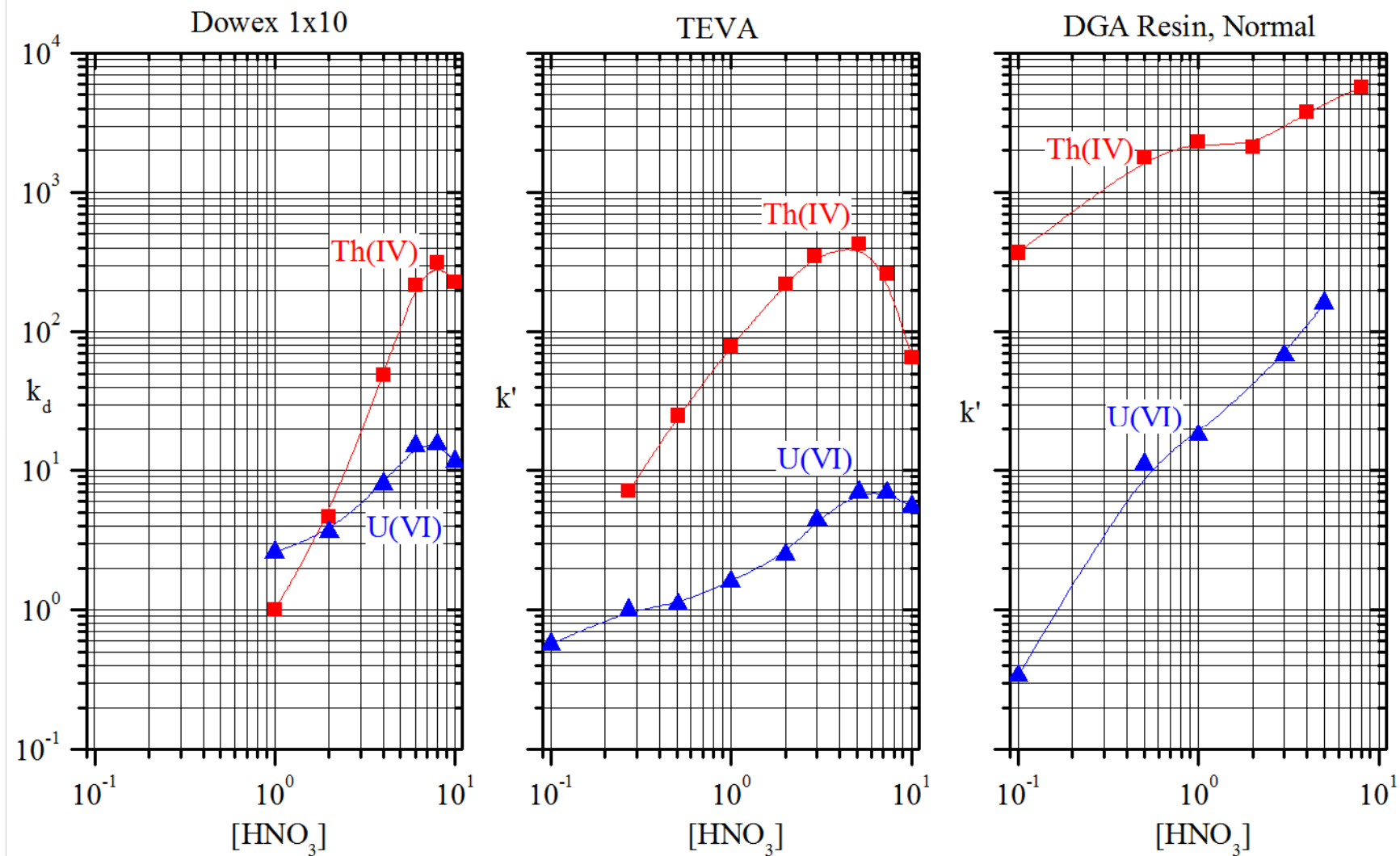

cpm Th-227 vs. Bed Volumes of Eluate

35 mL column of DGA Resin (12.6cm x 1.9cm), 100-150 μm , 18mL/min, 22(1) $^{\circ}\text{C}$

