

TANDEM SEPARATIONS OF RADIONUCLIDES USING AN AUTOMATED SYSTEM

Daniel R. McAlister and E. Philip Horwitz

55th Annual Radiobioassay and Radiochemical
Measurement Conference

October 26-30, 2009



Outline

- ARSIle Automated System


- Separation Challenges

- Examples of 2, 3, and 4 column systems

ARSIIe



Separation Challenges/Goals

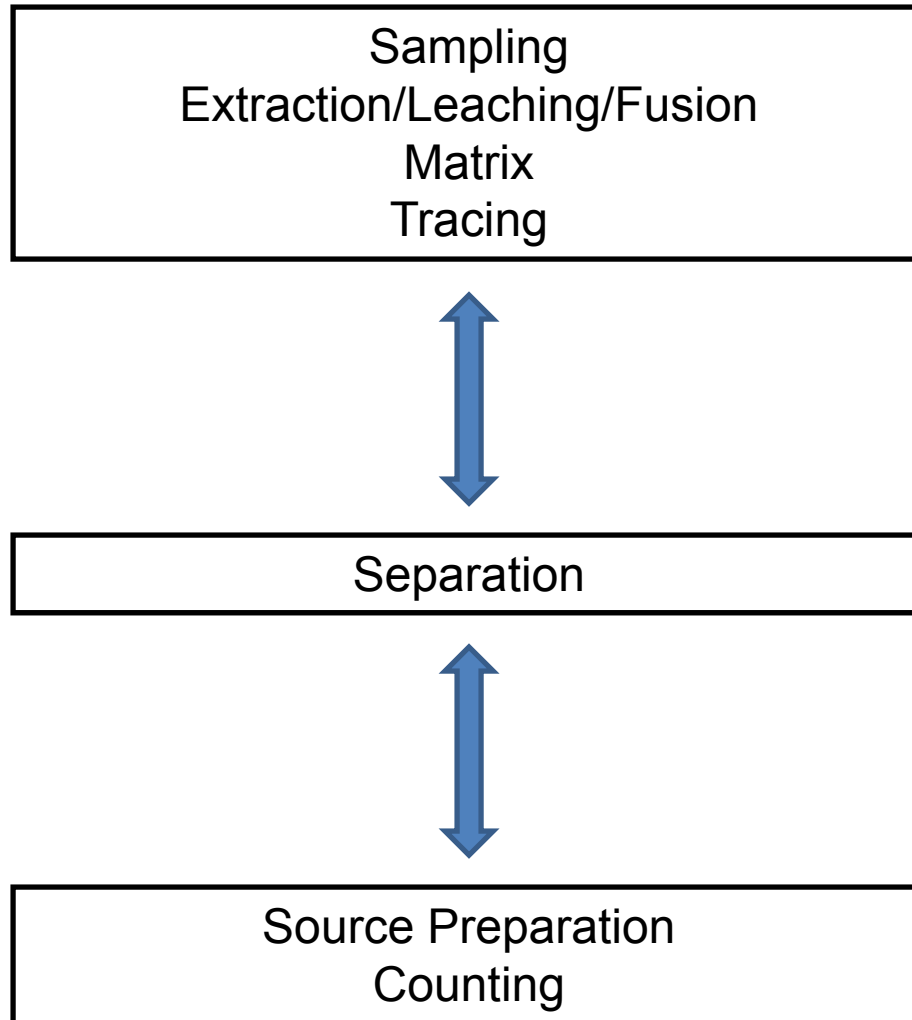

Separation Challenges/Goals

- Obtain desired analytes:
 - in the necessary purity
 - in a matrix which is compatible with the available measurement technique(s).
- Integrate system with sample extraction/fusion method
- Use as few different reagents as possible
- Use reagents that have long shelf-lives
- Use reagents that compatible with each other
- Minimize elution volumes

Available Tools

- TEVA

-Quaternary Amine: Th, Tc, Pu(IV), Np(IV), U(VI)

- UTEVA

-Phosphonate: Th, U(VI), Pu(IV), Np(IV), Po

- TRU

-Carbmoylphosphine oxide: Th, U(VI), Np(IV), Pu, Am

- DGA

-Diglycolamide: Th, U(VI), Np(IV), Pu, Am, ~Sr, ~Pb

- Sr Resin




-Crown Ether: Sr, Po, Pb, Pu(IV)

HNO₃

HCl

Both

Examples

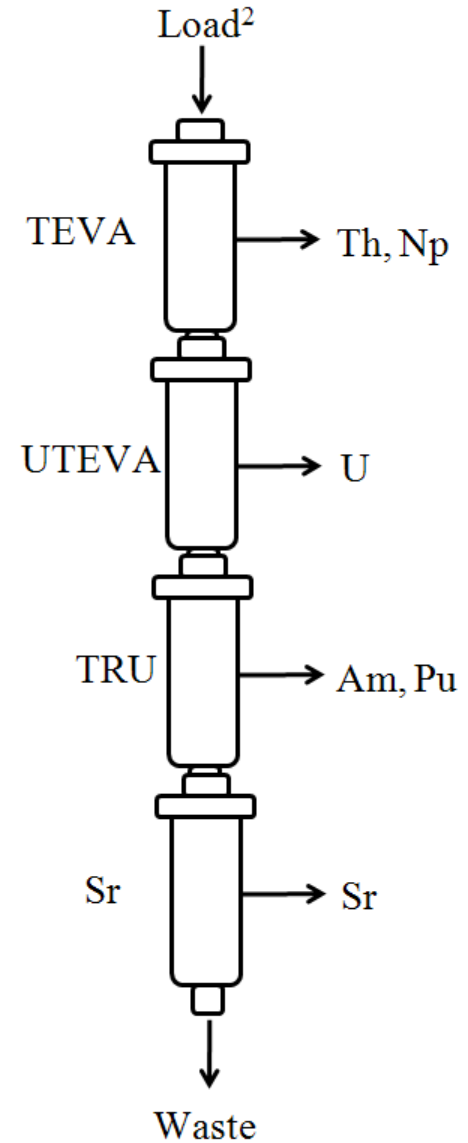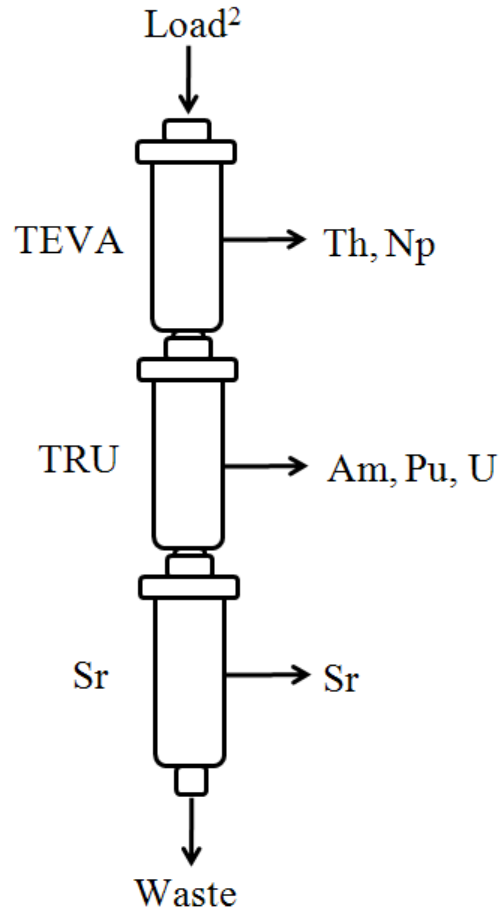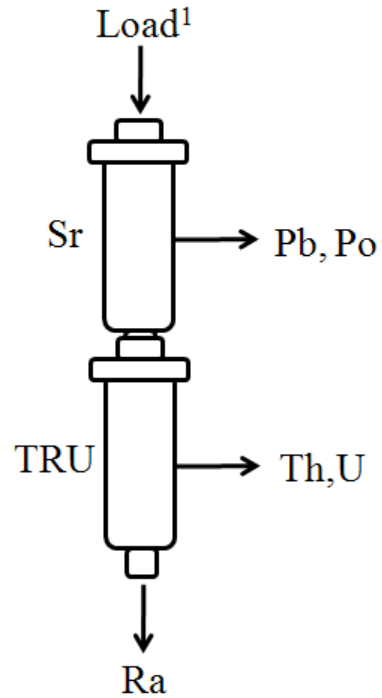


¹Load = 4M HCl + 0.1 M Ascorbic acid

²Load = 3M HNO₃ + 0.1M Ascorbic acid + 0.1 M Ferrous Sulfamate

1st Example: Pb, Po, U, Th, Ra

- NATO Project, Uranium Mining/Processing Sites in former Soviet Republics
- ^{238}U , ^{234}U , ^{230}Th , ^{226}Ra , ^{210}Pb and ^{210}Po from ^{238}U decay chain
- ^{232}Th , ^{228}Th , ^{224}Ra from ^{232}Th decay chain
- Soil samples leached with HCl
- Ascorbic acid added to limit interference from Fe(III)

Pb, Po, U, Th, Ra

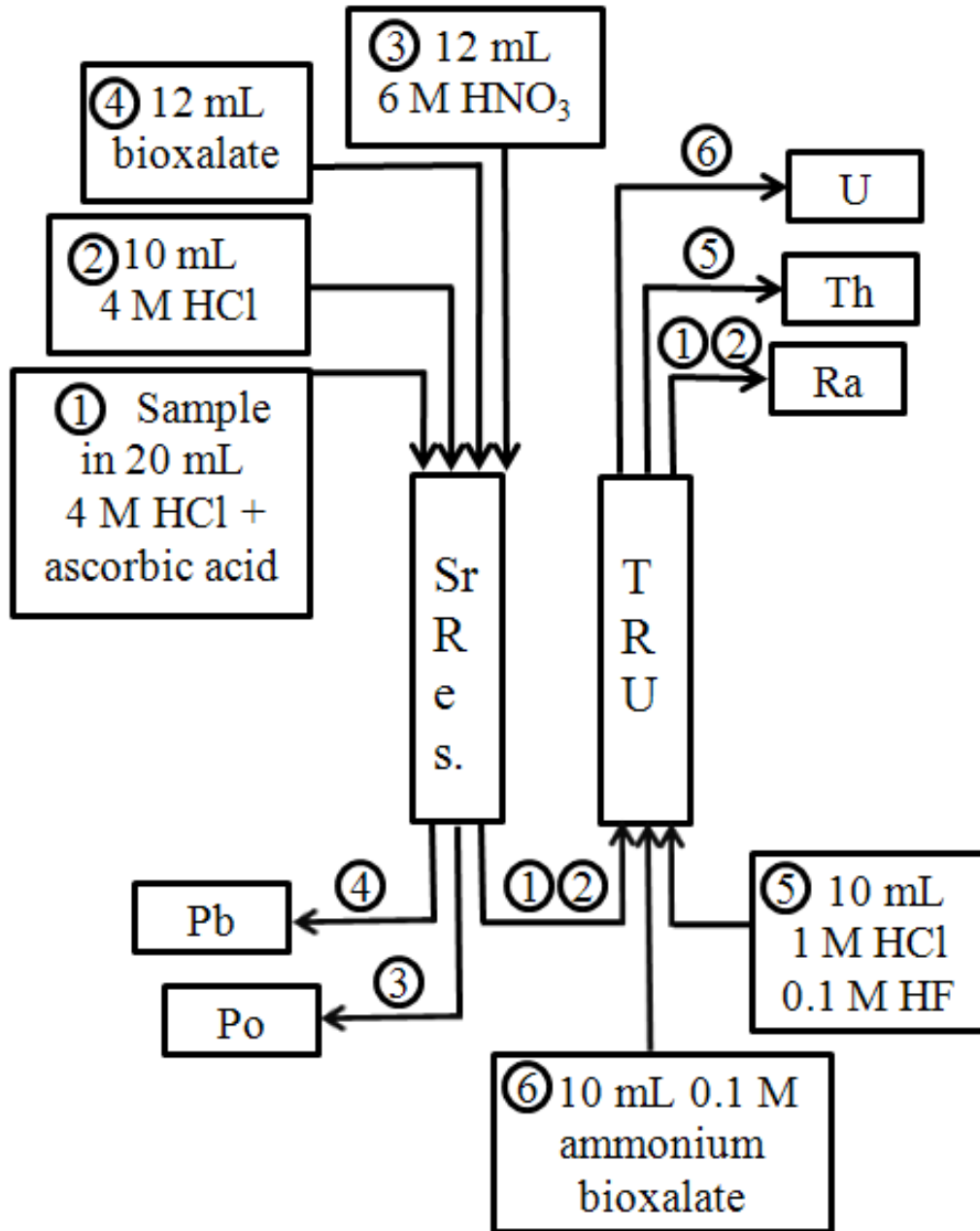

Direct precipitation or deposition of column eluate:

BaSO₄ precip → Ra

PbSO₄ precip → Pb


Autodeposition → Po

CeF₃ precip → U/Th

Followed by alpha spectroscopy


Pb, Po, U, Th, Ra

2.0 mL cartridge Sr Resin, 22(1)°C, 50-100 μm, 2 mL/min

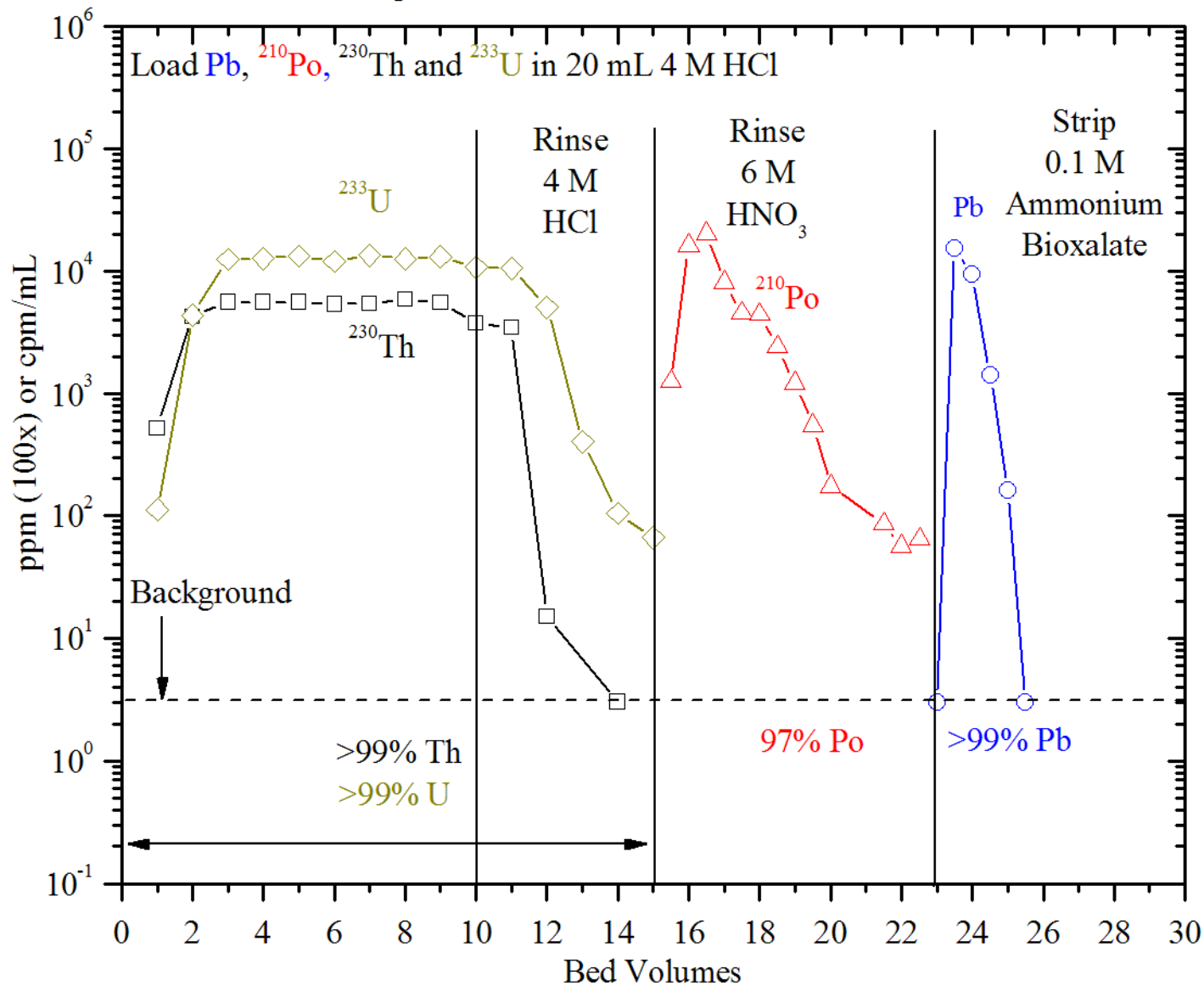


Pb, Po, U, Th, Ra


2.0 mL cartridge TRU, 22(1)°C, 50-100 μm, 2 mL/min

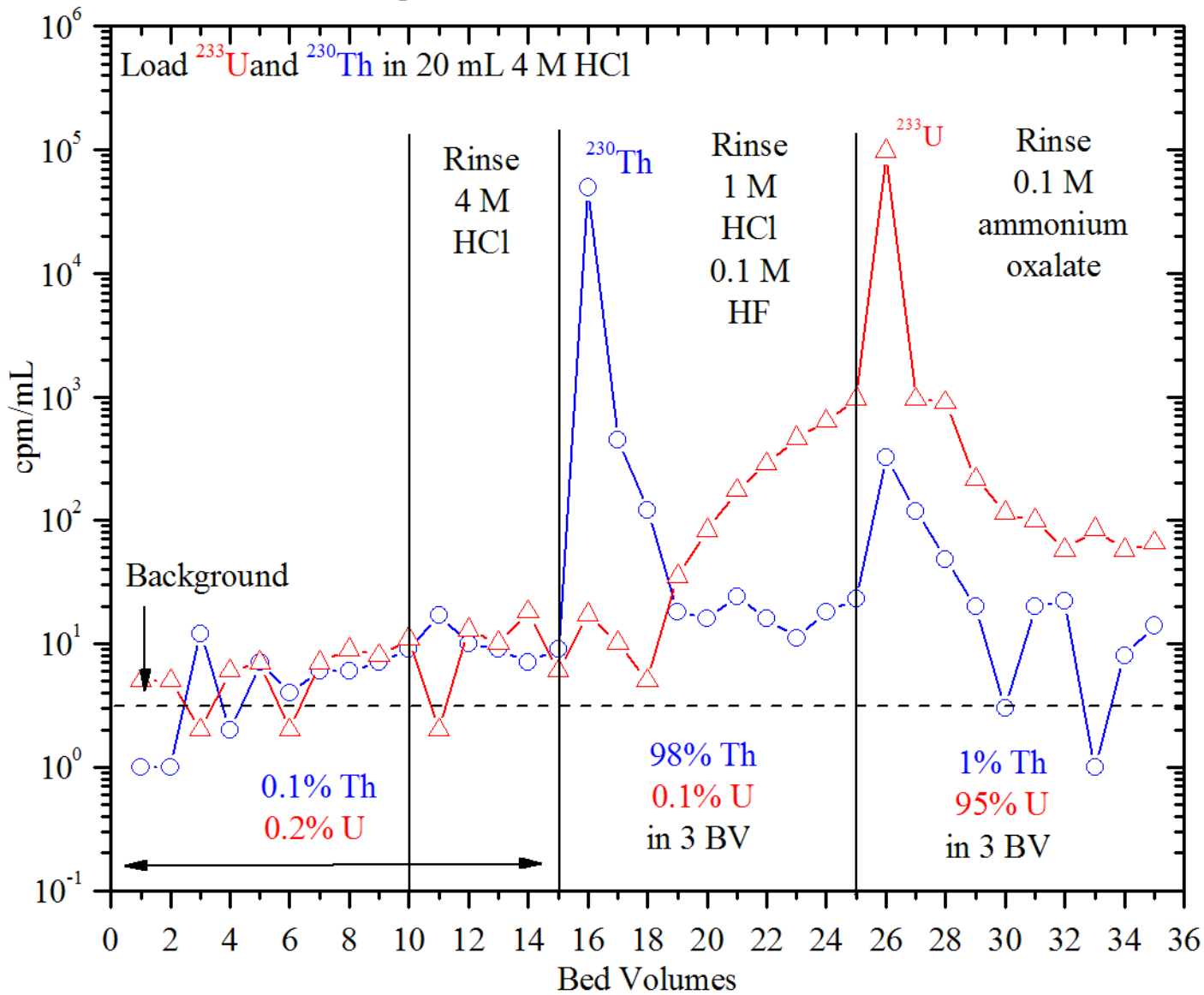


Pb, Po, U, Th, Ra

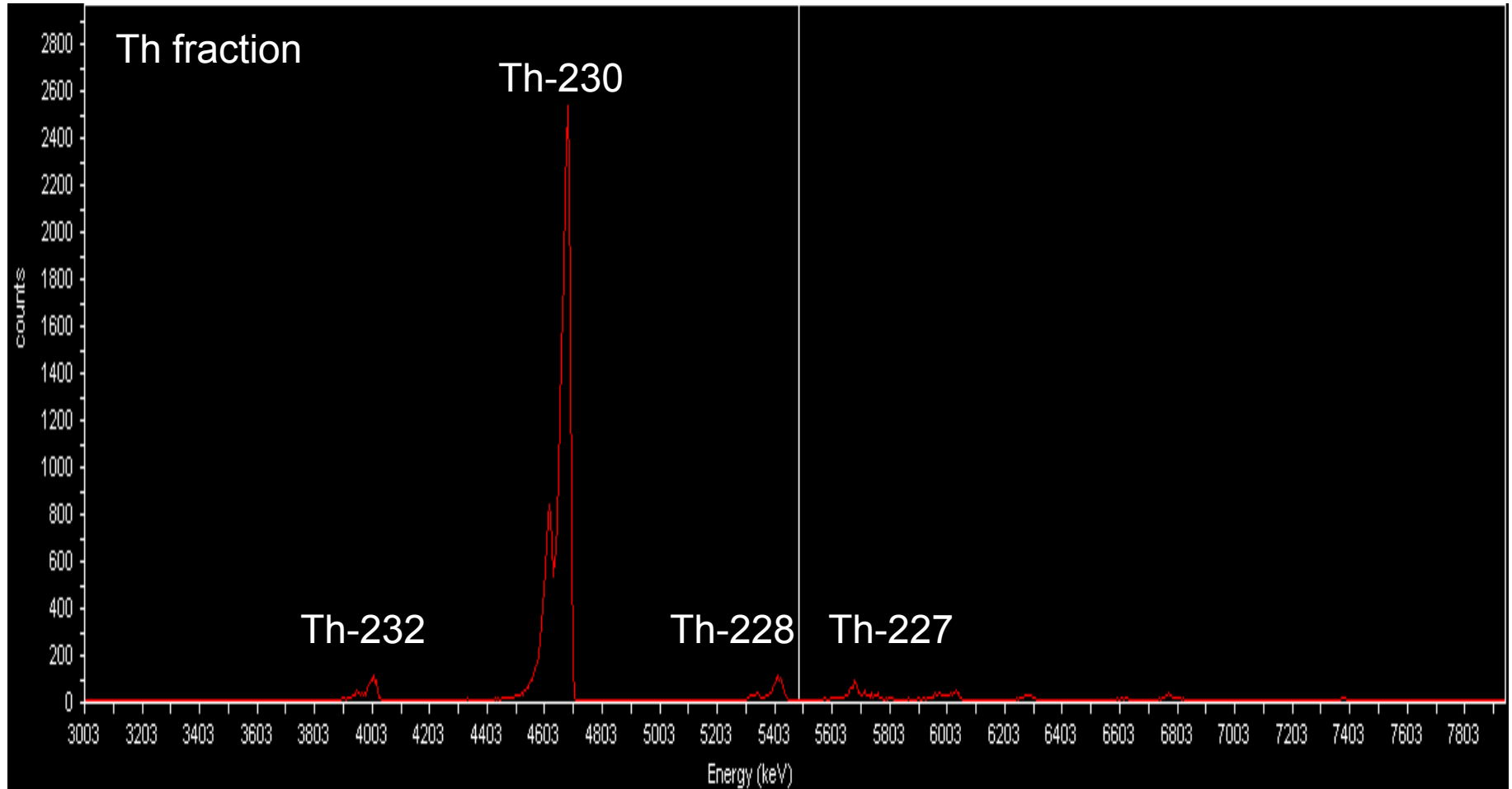


Pb, Po, U, Th, Ra

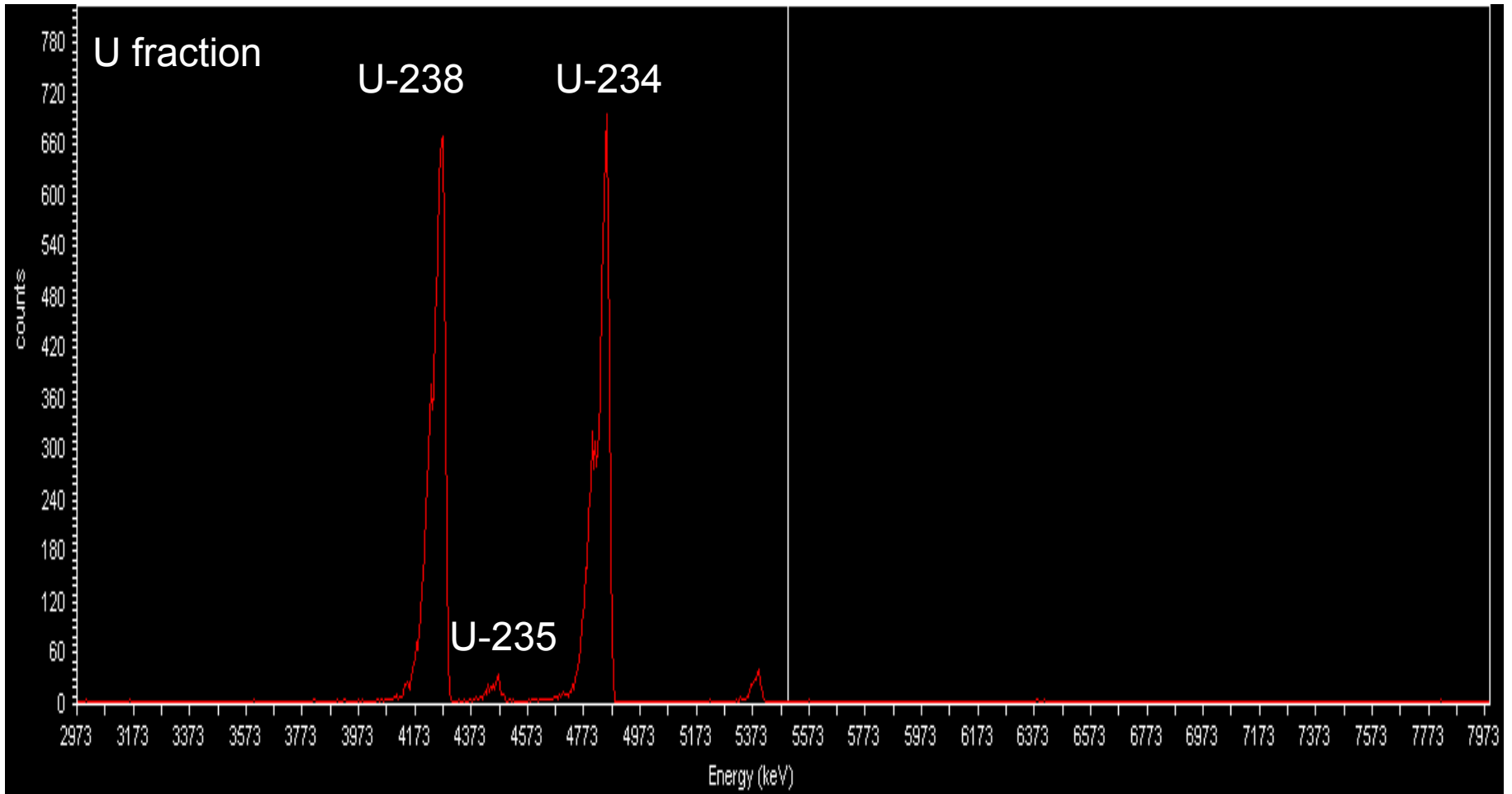

Pb, Po, U, Th, Ra

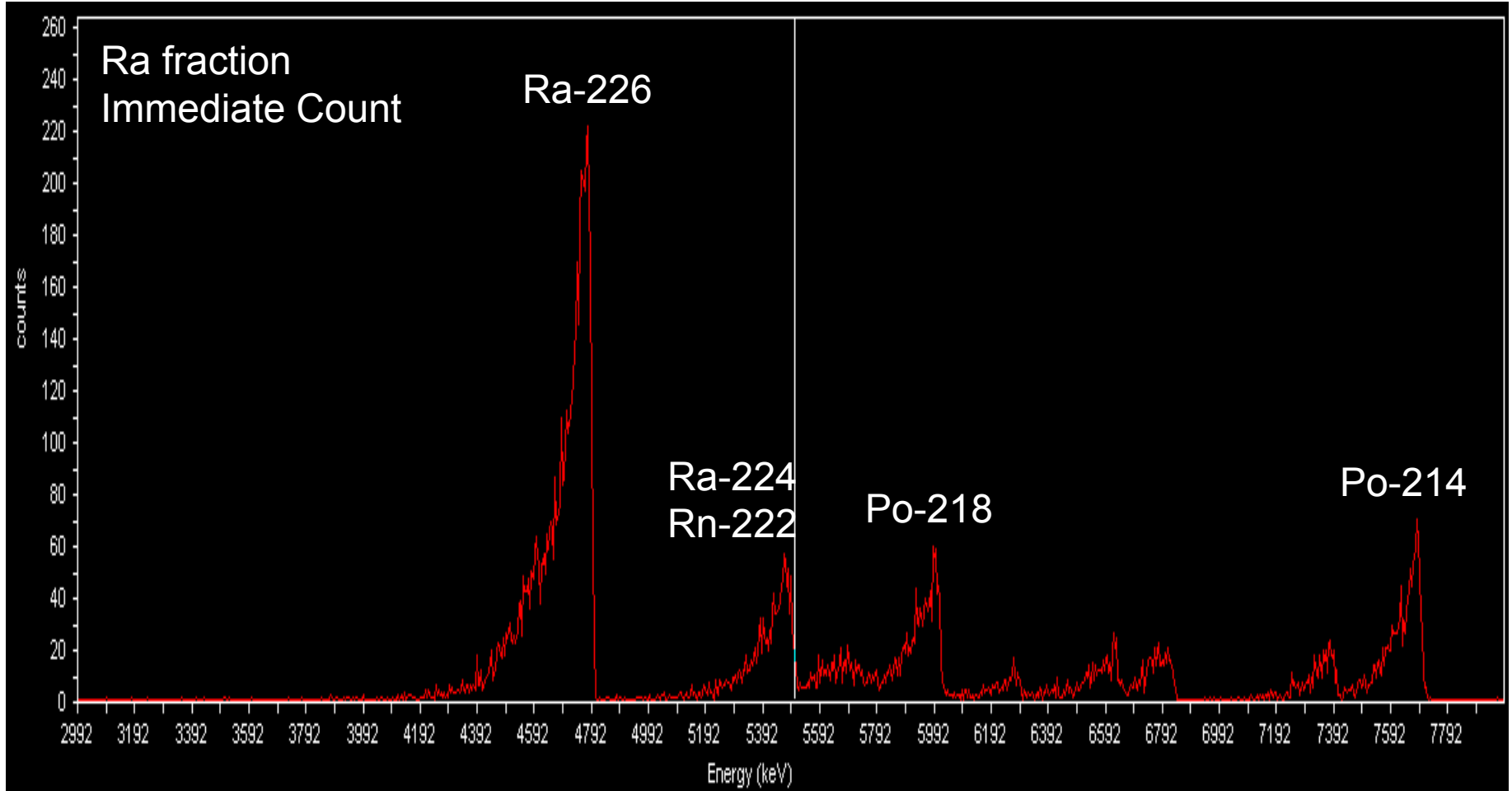


Pb, Po, U, Th, Ra

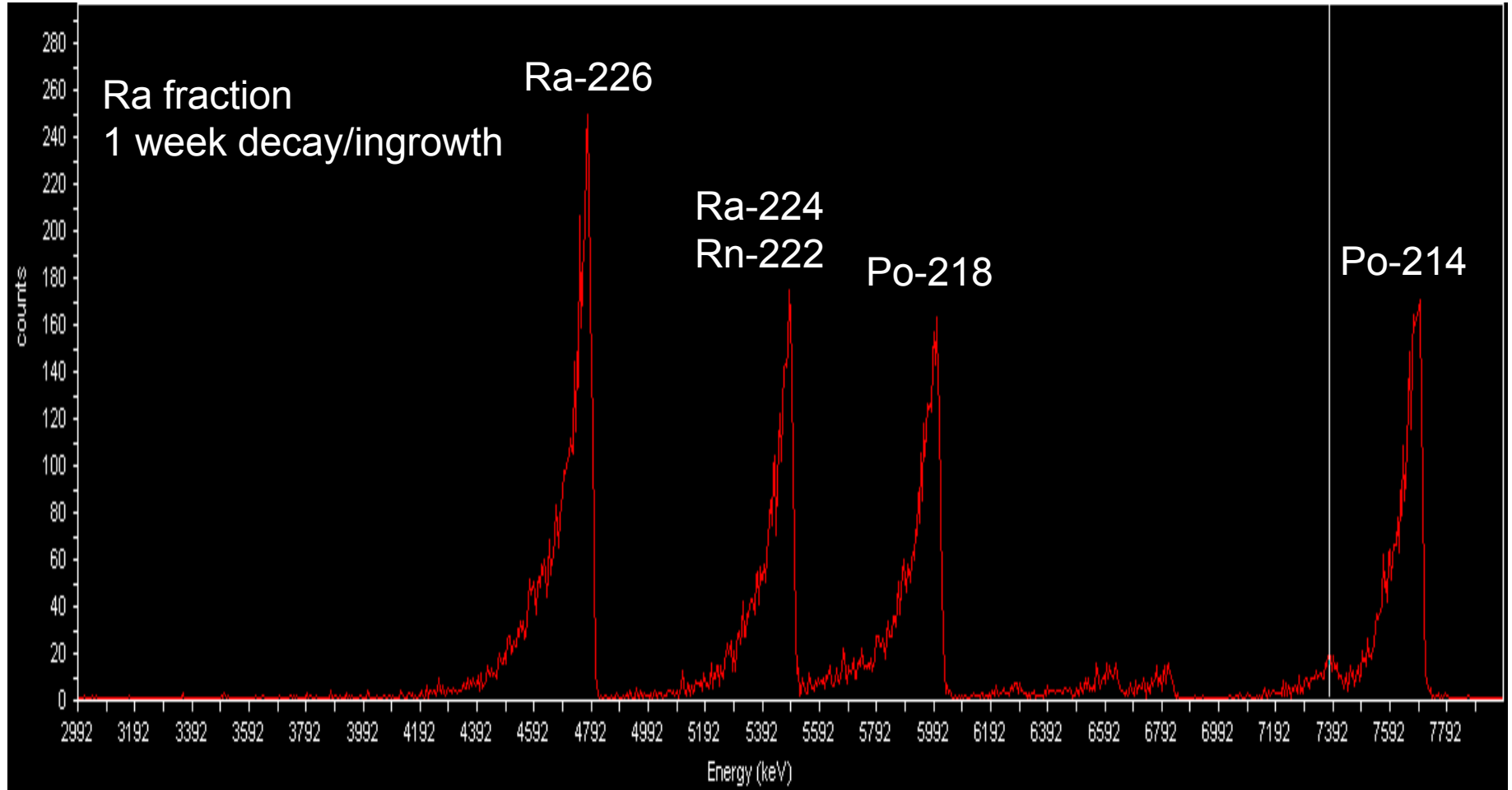

2nd Example: Th, Np, U, Am/Cm, Pu, Sr

- USAF Project (Marty Johnson Friday at 10:10)
- Soil samples, lithium borate fusion
- Dissolve Fusion with HNO_3
- Filter to remove silicates
- Adjust oxidation states with Ferrous Sulfamate/Ascorbic acid:
 - Np(IV)
 - Pu(III)
- Th, U, Am/Cm, Pu by alpha spectroscopy (CeF_3 precip.)
- Np by alpha, gas flow proportional and/or LSC (CeF_3 precip.)
- Sr by LSC or gas flow proportional

Th, Np, U, Am/Cm, Pu, Sr

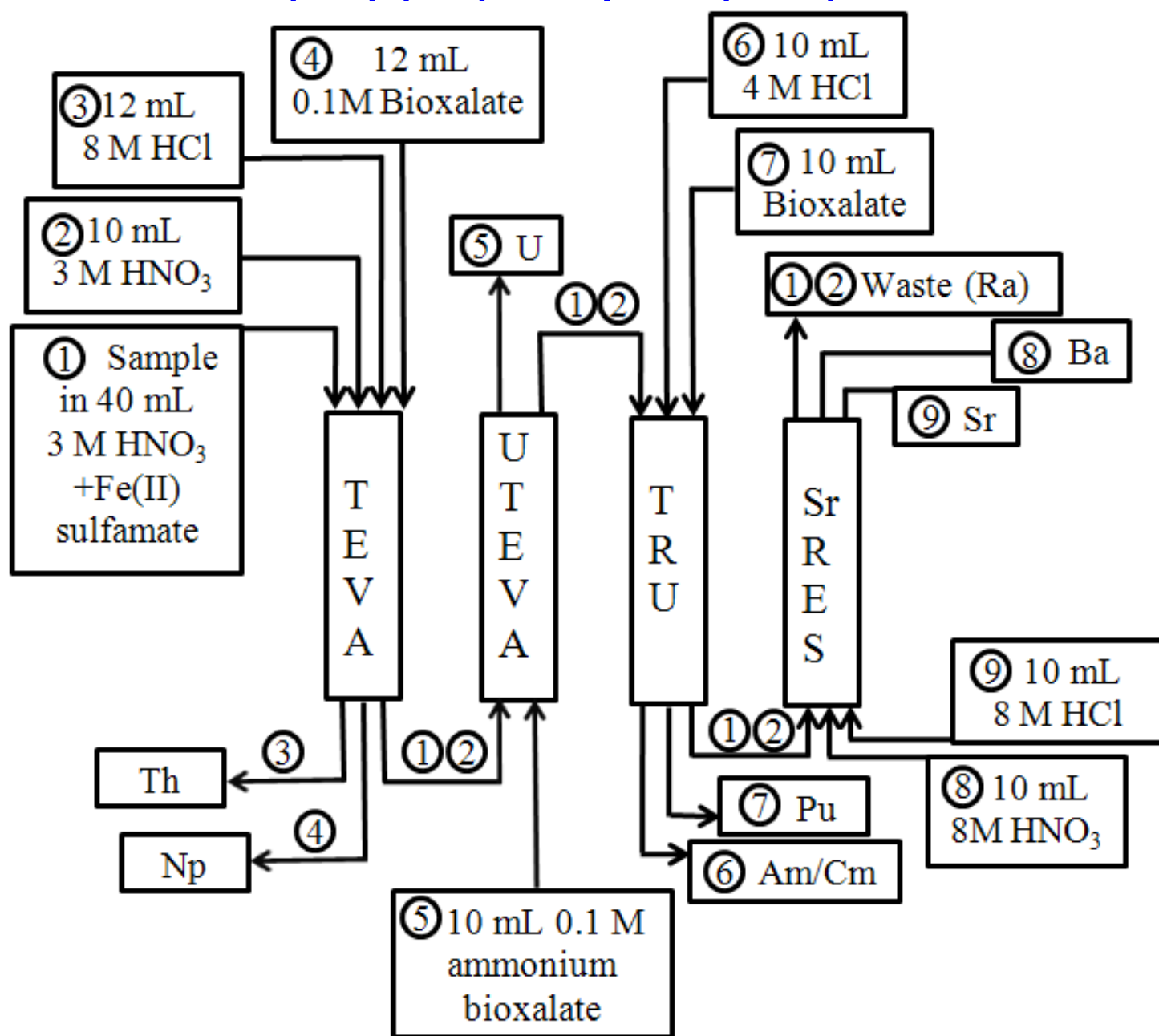

Th, Np, U, Am/Cm, Pu, Sr

Soil	Fraction	Nuclide	Tracer	
			Recovery (%) ^d	% Yield ^e
S2	TEVA-Th Strip	Th-228		
	12 mL	Th-229	219.7 ^f	
	8M HCl	Th-230		33.7 ^f
		Th-232		
S2	TEVA-Np 12 mL 0.1 M Biox	Np-237	N/A	
S2	UTEVA- U Strip	U-232	58.4	
	10 mL	U-233		
	0.1 M Bioxalate	U-234		98.6
		U-235		103.6
		U-238		109.8
S2	TRU-Am/Cm Strip	Am-241		123.5
	10 mL	Am-243	28.3	
	4M HCl	Cm-242		
		Cm-243		301.3
S2	TRU-Pu Strip	Pu-236		
	10 mL	Pu-238		127.8
	0.1 M Bioxalate	Pu-240		115.0
		Pu-242	62.7	
S2	Sr Resin Strip 10 mL 8 M HCl	Sr-90	86.8	

^aColumns: 2mL TEVA, UTEVA, TRU, Sr Resin (50-100 µm)

^bLoad: 45 mL 3M HNO₃ + Fe(II) sulfamate/ascorbic acid

^cRinse: 10 mL 3M HNO₃ + 0.1 M NaNO₂

^dAbsolute recovery of tracer through column chemistry and dissolution of melt

^eYield corrected for tracer recovery

^fPoor alpha spectrum resolution, overlap of Th-230 and Th-229 peaks

Th, Np, U, Am/Cm, Pu, Sr

Soil	Fraction	Nuclide	Tracer	
			Recovery (%) ^d	% Yield ^e
S3	TEVA-Th Strip	Th-228		
	12 mL	Th-229	86.2	
	8M HCl	Th-230		113.0
		Th-232		
S3	TEVA-Np 12 mL 0.1 M Biox	Np-237	N/A	
S3	UTEVA- U Strip	U-232	66.6	
	10 mL	U-233		
	0.1 M Bioxalate	U-234		102.7
		U-235		113.1
		U-238		102.2
S3	RE-2-Am/Cm Strip	Am-241		114.9
	10 mL	Am-243	55.9	
	4M HCl	Cm-242		
		Cm-243		94.0
S3	RE-2-Pu Strip	Pu-236		
	10 mL	Pu-238		87.2
	0.1 M Bioxalate	Pu-240		89.7
		Pu-242	97.6	
S3	Sr Resin Strip 10 mL 8 M HCl	Sr-90	107.0	

^aColumns: 2mL TEVA, UTEVA, RE-2, Sr Resin (50-100 µm)

^bLoad: 45 mL 3M HNO₃ + Fe(II) sulfamate/ascorbic acid

^cRinse: 10 mL 3M HNO₃ + 0.1 M NaNO₂

^dAbsolute recovery of tracer through column chemistry and dissolution of melt

^eYield corrected for tracer recovery

Conclusion

- Demonstrated two separation methods on the ARSIIe
 - 2 column (Po, Pb, U, Th, Ra)
 - 4 column (Th, Np, U, Am/Cm, Pu, Sr)
- Develop additional separations on ARSIIe
 - Single column methods (4 simultaneous samples)
 - Additional multiple column methods

Acknowledgment

Glenn Isensee, Mike Krummey, Lionel Johnson and Jim Harvey of Northstar Engineered Technologies for providing the ARSIIe instrument and training.

Marty Johnson for helpful discussions while developing the separation method.